


DOTS 2.0 – License activation


- Licenses are managed by admins
- Shipped as: .dotlicense file
- Must be activated to use the respective applications
 - Activate automatically if the DOTS instance has internet connection
 - Activate manually if the DOTS instance has no internet connection

Admin Page – Licenses tab

1. Log-in as admin
2. Click on the “A” in the upper right corner and select “Admin Settings

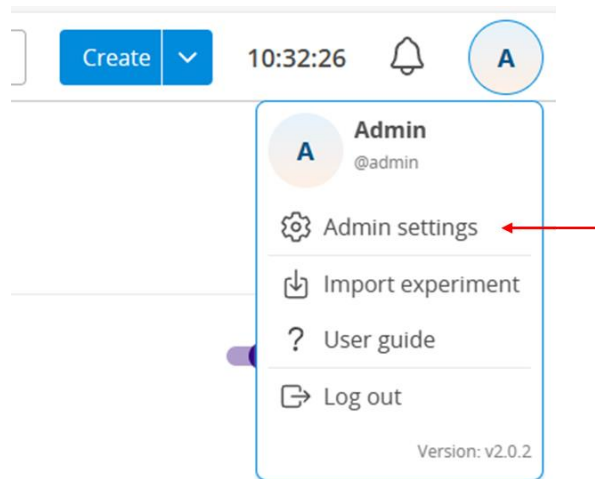


3. Click on the “Licenses” tab

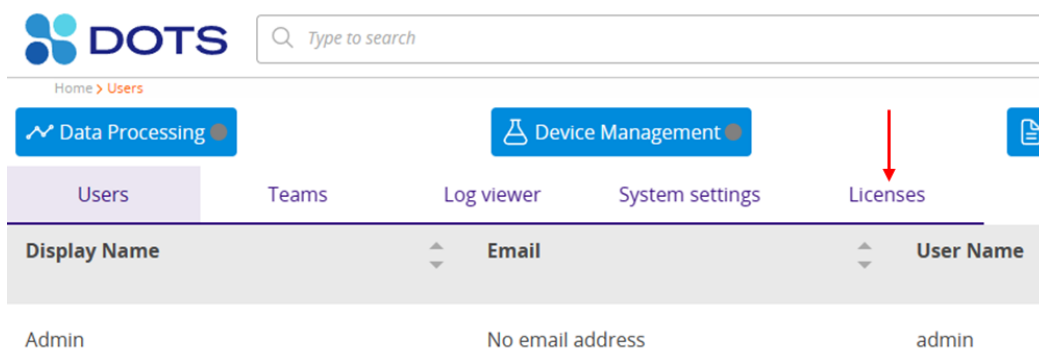


Online license activation (automatic)

1. Click on the blue “Add License” tab

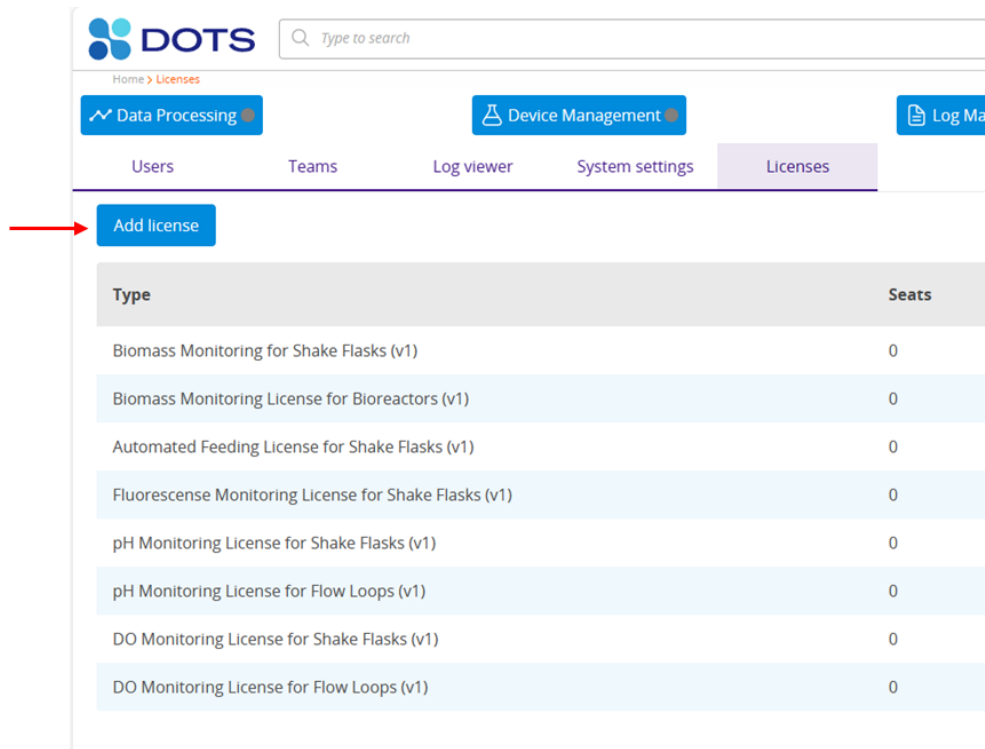

2. Upload your license file (that you got from sbi with your purchase). The seats contained in this license are shown
3. Click “Activate”. DOTS Software automatically connects to the Licensing Server and the licenses will get activated.

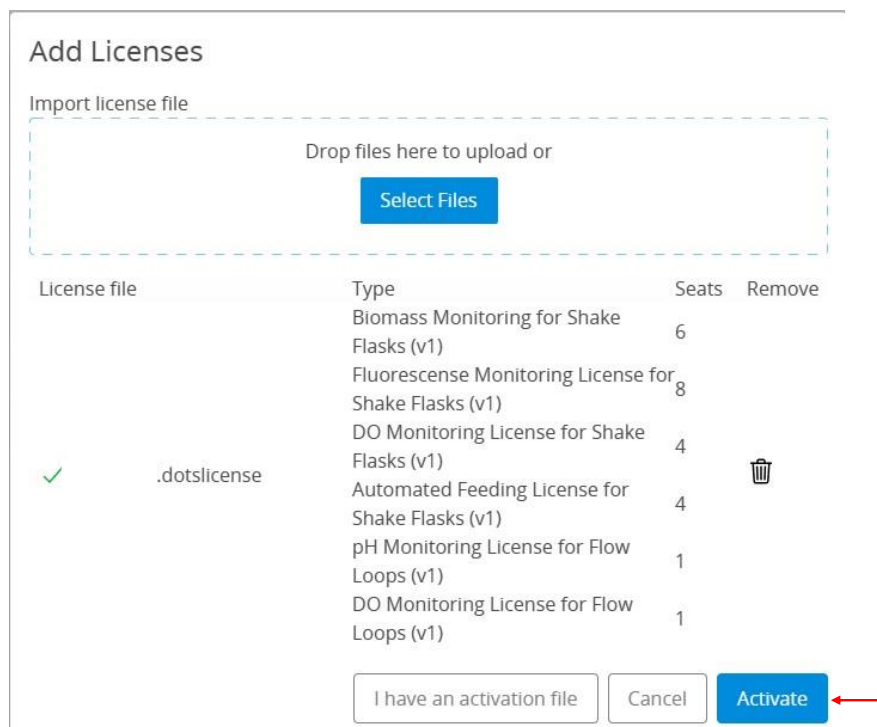

- Once the list of your activated Licenses is shown, click “Done”.

Add Licenses


License file	Type	Seats	Remove
	Biomass Monitoring for Shake Flasks (v1)	6	
	Fluorescence Monitoring License for Shake Flasks (v1)	8	
	DO Monitoring License for Shake Flasks (v1)	4	
✓ 2024-04-23T08_16_12.158Z.dots activation	Automated Feeding License for Shake Flasks (v1)	4	
	pH Monitoring License for Flow Loops (v1)	1	
	DO Monitoring License for Flow Loops (v1)	1	

Done

- You can now use the application templates you have the license for.

Offline license activation (manually)

- Click on the blue “Add License” tab

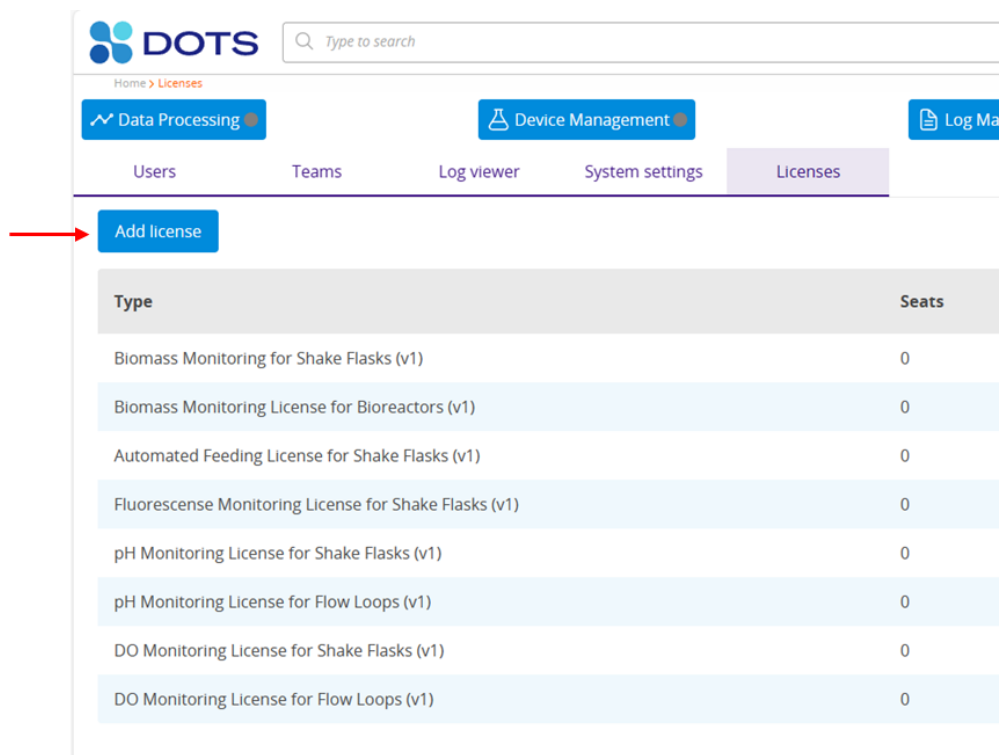

The screenshot shows the DOTS application interface. At the top, there is a search bar and navigation tabs: 'Data Processing', 'Device Management', and 'Log Ma'. Below these, there are sub-tabs: 'Users', 'Teams', 'Log viewer', 'System settings', and 'Licenses'. The 'Licenses' tab is selected. On the left side of the 'Licenses' tab, there is a blue 'Add license' button, which is highlighted by a red arrow. Below the button, there is a table with two columns: 'Type' and 'Seats'. The table lists various license types and their corresponding seat counts.

Type	Seats
Biomass Monitoring for Shake Flasks (v1)	0
Biomass Monitoring License for Bioreactors (v1)	0
Automated Feeding License for Shake Flasks (v1)	0
Fluorescence Monitoring License for Shake Flasks (v1)	0
pH Monitoring License for Shake Flasks (v1)	0
pH Monitoring License for Flow Loops (v1)	0
DO Monitoring License for Shake Flasks (v1)	0
DO Monitoring License for Flow Loops (v1)	0

2. DOTS activation server cannot be reached without internet connection. Click on “Download ID file”

Add Licenses

Activation server not reachable. Please download your machine's ID file and visit <https://licenses.scientificbio.com/activationKey/manual> to generate your activation file.

Once you have your activation key, press *I have my activation file*

Download ID file.

Cancel

I have my activation file

3. An automatic download of a .dotshwid file will be started. Transfer this file and the license (received from sbi) to a computer with internet connection.
4. Go to the website <https://licenses.scientificbio.com/activationKey/manual> to activate your license.

DOTS Activation

License files

Dateien auswählen
Keine Datei ausgewählt

This is the license file that you received from SBL.

Hardware ID file

Datei auswählen
Keine Datei ausgewählt

This file is generated by the DOTS Software when clicking "Download ID file".

Activate

Next steps

After clicking "Activate", the download of an Activation file ("*.dotsactivation") will be started. This file needs to be copied to a computer with access to the DOTS Software, for example via a USB drive. The Activation file can then be imported in the DOTS Software by opening the user interface, going to the Licenses tab in the Admin area, and following the steps after clicking the "Add License" button.

5. Upload your license file (received from sbi) and your hardware ID file (downloaded from the computer running DOTS). Click on “Activate”.

DOTS Activation

License files

Dateien auswählen
.dotslicense

This is the license file that you received from SBL.

Hardware ID file

Datei auswählen
DOTS-Unique-ID.dotshwid

This file is generated by the DOTS Software when clicking "Download ID file".

Activate

Next steps

After clicking "Activate", the download of an Activation file ("*.dotsactivation") will be started. This file needs to be copied to a computer with access to the DOTS Software, for example via a USB drive. The Activation file can then be imported in the DOTS Software by opening the user interface, going to the Licenses tab in the Admin area, and following the steps after clicking the "Add License" button.

6. An automatic download of a .dotsactivation file will start. Transfer this file to the computer that is running DOTS.
7. Repeat step 1 now clicking on “I have my activation file” in the popout.

Add Licenses

Activation server not reachable. Please download your machine's ID file and visit <https://licenses.scientificbio.com/activationKey/manual> to generate your activation file.

Once you have your activation key, press *I have my activation file*

Download ID file.

Cancel

I have my activation file

8. Upload the .dotsactivation file and click “Activate”.

Activate Licenses

Import activation file

Drop files here to upload or

Select Files

License file	Type	Seats	Remove
	Biomass Monitoring for Shake Flasks (v1)	6	
	Fluorescence Monitoring License for Shake Flasks (v1)	8	
	DO Monitoring License for Shake Flasks (v1)	4	
✓ 2024-04-23T08_16_12.158Z.dots activation	Automated Feeding License for Shake Flasks (v1)	4	
	pH Monitoring License for Flow Loops (v1)	1	
	DO Monitoring License for Flow Loops (v1)	1	

Cancel

Activate

9. Click “Done” to complete the activation of the license.
10. You can now use the application templates you have the license for.